



530 Main Street, Acton, MA 01720

Phone: (978)263-3584, **Fax:** (978)263-5086

Web Site: www.acton-research.com

Supplemental Instructions Indexable Slit / Aperture Assembly

Per customer specifications, this instrument is supplied with a special indexable slit aperture assembly. The slit assembly is bilaterally adjustable from 10 μ m to 3mm and operates as described in the SpectraPro[®] instruction manual. The aperture is 12mm high by 12mm wide. Either the slit or the aperture can be indexed into position under operating conditions by a control rod, located adjacent to the slit micrometer. When the control rod is up, the slit is in the operating position. When the control rod is down, the 12 x 12mm aperture is in the operating position.

NOTE: Slit width is not affected by movement of the control rod.